APPLICATIONS



Liquid-Liquid Extraction (LLE) vs. Novum[™] Simplified Liquid Extraction (SLE) Case Study: Analysis of Monomeric Phenolic Fermentation Inhibitors in Dilute-Acid Plant Hydrolysate

Ana B. Ibáñez and Stefan Bauer, Energy Biosciences Institute, University of California, Berkeley, CA 94720, USA

Introduction

Pretreatment of lignocellulosic biomass is a prerequisite in order to maximize glucose release from cellulose for subsequent fermentations or chemical conversions. A leading pretreatment involves dilute acid (e.g., 0.3 - 2 % sulfuric acid) at elevated temperatures (140 to 200 °C) with different holding times and biomass/acid loadings¹. A drawback of this procedure is the partial degradation of lignin, a polyphenolic macromolecule, to smaller phenolics that are released together with the sugars into the socalled hydrolysate. These phenolic compounds are toxic to microorganisms and are potential fermentation inhibitors ²⁻⁵.

Analyzing this phenolic profile is important when optimizing pretreatment or detoxification methods aiming to minimize inhibitor levels or to engineer more resistant microorganisms. In a widely employed sample preparation and cleanup step, the phenolic compounds undergo a liquid-liquid extraction using an organic solvent (e.g., ethyl acetate) and are then analyzed via GC/MS⁶. While effective, the liquid-liquid extraction can be time consuming, requiring 4 separate extractions and multiple centrifugation steps. In an effort to reduce time and sample handling steps, a simplified liquid extraction (SLE) using Novum SLE was employed, saving a significant amount of time and reducing the amount of required pipetting and vial handling steps.

Materials and Methods

Lignocellulosic biomass was prepared using two different methods, a traditional liquid-liquid extraction and simplified liquid extraction (SLE) using a Novum SLE MAX 96-well plate (part no. 8E-S138-5GA). Prior to each extraction method, the samples were pretreated as follows:

Pretreatment

A dilute-acid pretreatment hydrolysate (1.5 % sulfuric acid, 190 °C for 1 minute, 25 % biomass loading) from *Miscanthus X giganteus* was filtered using a PES (polyethersulfone) 0.45 μ m filter and was spiked with 20 μ g/mL iso-propylphenol as an internal standard (IS). The pretreated samples were then extracted by either liquid-liquid extraction or SLE (below).

Liquid-Liquid Extraction Procedure

1 mL aliquots of the spiked hydrolysate from the pretreatment step were extracted 4x each with 0.5 mL of ethyl acetate. After each extraction, a centrifugation step (1500 g for 1 minute) was applied for phase separation. The combined extracts of each sample were dried over sodium sulfate⁶ (samples A).

Simplified Liquid Extraction (SLE) Procedure

 $400\,\mu$ L aliquots of the spiked hydrolysate from the pretreatment step were loaded into the wells of a Novum SLE MAX 96-well plate. The sample was allowed to soak into the sorbent for 5 minutes and then the ethyl acetate was applied to elute the analytes of interest. Three different volumes were applied to determine the most effective elution volume. Either $1 \times 0.4 \text{ mL}$ (samples B), $1 \times 0.8 \text{ mL}$ (samples C), or $2 \times 0.4 \text{ mL}$ (samples D) of ethyl acetate was applied using a short pulse of vacuum. No further drying of ethyl acetate was performed after the elution.

GC/MS Conditions

Prior to GC analysis, 100 µL aliquots of the extracts (samples A -D) were incubated with 50 µL BSTFA at 70 °C for 30 minutes and then analyzed by GC/MS⁶. All extractions were performed in triplicate and replicates (1-3) were run in the following order: samples A1 - D1, then A2 - D2, then A3 - D3 to avoid any bias because of potential changes in sample composition of unprocessed samples due to the long run time (67 minutes). 1 µL was injected in splitless mode onto a 5-ms capillary GC column (30m x 0.25 mm x 0.25 µm). An Agilent[®] 7890A gas chromatograph coupled to an Agilent 5975C single quadrupole mass spectrometer was used for analysis with the following settings: Injector and transfer line temperature 280 °C, carrier gas: helium at 1 mL/min, temperature program: 75 °C for 3 min isocratic, 5 °C/min to 150 °C, 0.5 °C/min to 160°C, 2°C/min to 190°C, 5°C/min to 240°C, 70°C/min to 325 °C. 3 min isocratic. Ions were detected in full scan mode m/z 35-500.

Results and Discussion

Figure 1 shows the relative recoveries of 25 selected phenolic compounds from hydrolysate eluted from the Novum SLE MAX 96-well plate using the 3 elution procedures (0.4 mL, 0.8 mL and 2x 0.4 mL) compared to the values obtained by the conventional liquid-liquid extraction (=100% on the representative chart). Overall the Novum SLE extraction produced recoveries that were comparable to the traditional liquid-liquid extraction procedure in a fraction of the time (a five-fold decrease in the amount of sample preparation time). The 0.4 mL ethyl acetate elution resulted in a general lower recovery of the analytes compared to the 0.8 mL and 2x 0.4 mL volume extractions. Exceptions were 4-hydroxy-3-methoxy-cinnamaldehyde (Figure 1, analyte 9) with almost identical recovery levels observed (95% and 93%, respectively) for all elution volumes and sinapaldehyde (Figure 1, analyte 12) with a higher recovery for the 0.4 mL extraction volume (93 %) compared to the other extraction volumes applied (89 % and 85 %, respectively). However, a 0.4 mL extraction volume was generally deemed to be suboptimal. Increasing the elution volume to 0.8 mL improved the recovery with most recoveries reaching >90%. Slightly lower yields were observed for 4-hydroxybenzaldehyde (87 %) (Figure 1, analyte 1) and vanillin (89 %) (Figure 1, analyte 2). Only the homologs 2-guaiacylacetaldehyde (Figure 1, analyte 14) and 2-syringacetaldehyde (Figure 1, analyte 16) exhibited lower recovery levels of 75-76% and 73-75%, respectively. Applying the 0.8 mL elution volume in two steps (2x 0.4 mL) did not improve the recovery as compared to a single aliquot of 0.8 mL.

APPLICATIONS



Conclusion

When extracting 0.4 mL of a dilute-acid plant hydrolysate using a Novum[™] SLE MAX 96-well plate, a 0.8 mL elution volume of ethyl acetate resulted in high recoveries of >90 % for most phenolic analytes. No further drying of the ethyl acetate phase was necessary to ensure successful derivatization for subsequent GC/MS analysis. Compared to the conventional liquid-liquid extraction, we estimate an up to five-fold decrease in sample preparation time. Furthermore, the Novum SLE extraction also greatly simplifies sample preparation by reducing the amount of required pipetting and vial handling steps. Therefore, sample preparation using the Novum SLE resulted in a phenolic inhibitor profile that was very similar and comparable to the conventional liquid-liquid extraction but at a fraction of the time and labor.

Figure 1. Relative recovery (%) after extraction using Novum SLE under various extraction solvent volumes as compared to a liquid-liquid extraction

References

- Humbird D, Davis R, Tao L, Kinchin C, Hsu D, Aden A, Schoen P, Lukas J, Olthof B, Worley M, Sexton D, Dudgeon D (2011) Process Design and Economics for Biochemical Conversion of Lignocellulosic Biomass to Ethanol: Dilute-Acid Pretreatment and Enzymatic Hydrolysis of Corn Stover. Technical Report NREL/TP-5100-47764. United States National Renewable Energy Laboratory, US Department of Energy. May 2011
- Jönsson LJ, Palmqvist E, Nilvebrant N-O, Hahn-Hägerdal B (1998) Detoxification of wood hydrolysates with laccase and peroxidase from the white-rot fungus Trametes versicolor. Appl Microbiol Biotechnol 49:691-697
- Luo C, Brink DL, Blanch HW (2002) Identification of potential fermentation inhibitors in conversion of hybrid poplar hydrolyzate to ethanol. Biomass Bioenergy 22:125-138
- Larsson S, Reimann A, Nilvebrant N-O, Jönsson LJ (1999) Comparison of different methods for the detoxification of lignocellulose hydrolyzates of spruce. App Biochem Biotechnol 77-79:91- 103
- Buchert J, Niemelä K, Puls J, Poutanen K (1990) Improvement in the fermentability of steamed hemicellulose hydrolysate by ion exclusion. Process Biochem Int 176-180
- Mitchell VD, Taylor CM, Bauer S (2014) Comprehensive analysis of monomeric phenolics in dilute acid plant hydrolysates. Bioenerg Res 7:654-669



Analyte Key:		
1	4-hydroxybenzaldehyde	
2	vanillin	
3	acetovanillone	
4	4-hydroxybenzoic acid	
5	syringaldehyde	
6	vanillic acid	
7	homovanillic acid	
8	3,4-dihydroxybenzoic acid	
9	4-hydroxy-3-methoxy-cinnamaldehyde	
10	syringic acid	
11	trans-p-coumaric acid	
12	sinapaldehyde	

Analyte Key cont'd:

13	trans-ferulic acid	
14	2-guaiacylacetaldehyde	
15	guaiacylacetone	
16	2-syringacetaldehyde	
17	3-guaiacylpropanol	
18	2-hydroxy-1-guaiacylpropanone	
19	3-guaiacylacetol	
20	3-syringylacetol	
21	1-guaiacylacetol	
22	3-(4-hydroxyphenyl)acetol	
23	1-syringylacetol	
24	2-hydroxy-1-guaiacylethanone	
25	2-hydroxy-1-syringylethanone	



PLICATIONS

Ordering Information

Novum[™] Simplified Liquid Extraction (SLE) 96-Well Plates Description Unit Part No. 8E-S138-FGA Novum SLE MINI 96-Well Plate 1/pk 8E-S138-5GA Novum SLE MAX 96-Well Plate



96-Well Plate Accessories				
Part No.	Description	Unit		
Collection Plates (deep well, polypropylene)				
AH0-7192	96-Well Collection Plate, 350 µL/well	50/pk		
AH0-7193	96-Well Collection Plate, 1 mL/well	50/pk		
AH0-7194	96-Well Collection Plate, 2 mL/well	50/pk		
AH0-8635	96-Well Collection Plate, 2 mL/well Square/Round-Conical	50/pk		
AH0-8636	96-Well Collection Plate, 2 mL/well Round/Round, 8 mm	50/pk		
AH0-7279	96-Well Collection Plate, 1 mL/well Round, 7 mm	50/pk		
Sealing Mats				
AH0-8597	Sealing Mats, Pierceable, 96-Square Well, Silicone	50/pk		
AH0-8598	Sealing Mats, Pre-Slit, 96-Square Well, Silicone	50/pk		
AH0-8631	Sealing Mats, Pierceable, 96-Round Well 7 mm, Silicone	50/pk		
AH0-8632	Sealing Mats, Pre-Slit, 96-Round Well 7 mm, Silicone	50/pk		
AH0-8633	Sealing Mats, Pierceable, 96-Round Well 8 mm, Silicone	50/pk		
AH0-8634	Sealing Mats, Pre-Slit, 96-Round Well 8 mm, Silicone	50/pk		
AH0-7362	Sealing Tape Pad	10/pk		
Vacuum Manifold				
AH0-8950	96-Well Plate Manifold, Universal with Vacuum Gauge	ea		

1/pk

Novum SLE Tubes

Part No.	Description	Unit
8B-S138-FAK	Novum SLE 1 cc tubes	100/pk
8B-S138-5BJ	Novum SLE 3 cc tubes	50/pk
8B-S138-JCH	Novum SLE 6 cc tubes	30/pk
8B-S138-KDG	Novum SLE 12 cc tubes	20/pk

Tube Accessories

Vacuum Manifolds			
Part No.	Description	Unit	
AH0-6023	12-Position Vacuum Manifold Set	ea	
AH0-6024	24-Position Vacuum Manifold Set	ea	



If Novum SLE products do not perform as well or better than your current SLE product, return the product with comparative data within 45 days for a FULL REFUND.

PI ICATIONS



Australia

t: +61 (0)2-9428-6444 f: +61 (0)2-9428-6445 auinfo@phenomenex.com

Austria

- t: +43 (0)1-319-1301 f: +43 (0)1-319-1300
- anfrage@phenomenex.com

Belaium

t: +32 (0)2 503 4015 (French) t: +32 (0)2 511 8666 (Dutch) f: +31 (0)30-2383749 beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681 f: +1 (310) 328-7768

info@phenomenex.com

China t: +86 (0)20 2282-6668

f: +86 (0)20 2809-8130

chinainfo@phenomenex.com

Denmark t: +45 4824 8048

f: +45 4810 6265 nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063 f: +45 4810 6265 nordicinfo@phenomenex.com

France t: +33 (0)1 30 09 21 10 f: +33 (0)1 30 09 21 11

franceinfo@phenomenex.com

- **Germany** t: +49 (0)6021-58830-0 f: +49 (0)6021-58830-11
- anfrage@phenomenex.com

India

t: +91 (0)40-3012 2400 f: +91 (0)40-3012 2411 indiainfo@phenomenex.com

Ireland

- t: +353 (0)1 247 5405
- f: +44 1625-501796 eireinfo@phenomenex.com

Italv

- t: +39 051 6327511 f: +39 051 6327555
- italiainfo@phenomenex.com

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department at international@phenomenex.com

Luxembourg t: +31 (0)30-2418700

f: +31 (0)30-2383749 nlinfo@phenomenex.com

Mexico t: 001-800-844-5226

- f: 001-310-328-7768
- tecnicomx@phenomenex.com

The Netherlands t: +31 (0)30-2418700 f: +31 (0)30-2383749

- nlinfo@phenomenex.com

New Zealand t: +64 (0)9-4780951

- f: +64 (0)9-4780952
- nzinfo@phenomenex.com

Norway t: +47 810 02 005

- f: +45 4810 6265
- nordicinfo@phenomenex.com

- Puerto Rico t: +1 (800) 541-HPLC f +1 (310) 328-7768
- info@phenomenex.com

Spain

- t: +34 91-413-8613 f: +34 91-413-2290 espinfo@phenomenex.com

Sweden t: +46 (0)8 611 6950

- f: +45 4810 6265
- nordicinfo@phenomenex.com

United Kingdom

- t: +44 (0)1625-501367 f: +44 (0)1625-501796
- ukinfo@phenomenex.com

USA

- t: +1 (310) 212-0555
- f: +1 (310) 328-7768 info@phenomenex.com

All other countries Corporate Office USA t: +1 (310) 212-0555

- f: +1 (310) 328-7768
- info@phenomenex.com

Terms and Conditions

Subject to Phenomenex Standard Terms and Conditions which may be viewed at www.phenomenex.com/TermsAndConditions.

Trademarks

Novum is a trademark of Phenomenex. Agilent is a registered trademark of Agilent Technologies. Disclaimer

UC Berkeley, including its Energy Biosciences Institute (EBI), is not affiliated with Phenomenex and does not support, endorse, or sponsor it or its products/services.

Novum is patent pending.

© 2015 Phenomenex, Inc. All rights reserved.

