
Phenomenex ...breaking with tradition

APPLICATIONS

Related Substances Assay of the Drug Product Clarithromycin on a Kinetex 3.5 µm XB-C18 Column According to Ph. Eur. Monograph 1651

Matthew Trass¹, Simon Lomas¹ and Heiko Behr²

¹Phenomenex, Inc., 411 Madrid Ave., Torrance, CA 90501 USA
²Phenomenex, Ltd., Zeppelinstr. 5, 63741 Aschaffenburg, Germany

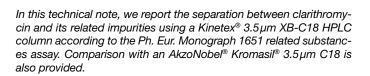

Introduction

The European Pharmacopoeia (Pharmacopoea Europaea, Ph. Eur.) is a single reference work for the quality control of medicines in the signatory states of the Convention on its elaboration. The official standards published within provide a legal and scientific basis for quality control during the development, production, and marketing processes. They concern the qualitative and quantitative composition and the tests to be carried out on medicines, on the raw materials used in production of medicines and on the intermediates of synthesis. All producers of medicines and/or substances for pharmaceutical use must therefore apply these quality standards in order to market their products.

Clarithromycin is an antibiotic used to treat a variety of bacterial infections including strep throat, pneumonia, skin infections and Lyme disease.

Figure 1. Clarithromycin



Dr. Heiko Behr

European Pharmaceutical Senior Business Development Manager

In addition to chromatography, Dr. Behr has an absolute love for heavy metal music and skiing in the Alps.







Materials and Methods

The analysis was performed using an Agilent® 1100 equipped with a multiple wavelength length UV detector. The Kinetex 3.5 µm XB-C18 column used for analysis was obtained from Phenomenex (Torrance, CA, USA). The Kromasil 3.5 µm C18 used for analysis was obtained from AkzoNobel, Separation Products (SE-445 80 Bohus, Sweden). All solvents were purchased from Honeywell (Morristown, NJ, USA) and Sigma-Aldrich (St. Louis, MO, USA). Ph. Eur. Standard Clarithromycin for peak identification CRS Y0000321 was purchased from European Directorate for the Quality of Medicines & HealthCare (EDQM) − Council of Europe; Postal address: 7 Allée Kastner CS 30026 F - 67081 STRASBOURG (France).

Ph. Eur. Monograph 1651

Ph. Eur. Monograph 1651 system suitability reference solution D. Dissolve 15 mg of Clarithromycin for peak identification CRS in 5 mL of acetonitrile and dilute to 10 mL with water.

Column:

- size: I = 0.10 m, $\emptyset = 4.6 mm$
- stationary phase: octadecylsilyl silica gel for chromatography R (3.5 μm)
- temperature: 40°C

Mobile phase:

- A: a 4.76 g/L solution of potassium dihydrogen phosphate adjusted to pH 4.4 with dilute phosphoric acid
- B: Acetonitrile

 $\textit{Gradient:}\ 25\,\%$ mobile phase B to 60 % mobile phase B in 32 min, hold for 2 min

Flow rate: 1.1 mL/min

Detection: spectrophotometer at 205 nm

Injection: 10 µL

Relative retention with reference to Clarithromycin (retention time = about 11 min): impurity A = about 0.42; impurity J = about 0.63; impurity L = about 0.74; impurity L = ab

System suitability: reference solution D:

 peak-to-valley ratio: minimum 3.0, where Hp= height above the baseline of the peak due to impurity D and Hv = height above the baseline of the lowest point of the curve separating this peak from the peak due to clarithromycin in the chromatogram obtained with reference solution D.



Result and Discussion

In this tech note we report the separation between Clarithromycin (the active pharmaceutical ingredient) and its related impurities per the related substances section of Ph. Eur. Monograph 1651, as shown in **Figure 2**. On the Kinetex column the Clarithromycin peak elutes at 9.25 min. The peak-to-valley ratio between impurity D and clarithromycin is 10.8 which passes the system suitability minimum requirement of 3.0. As a comparison, a Kromasil® 3.5 µm C18 was also used for the analysis, as shown in **Figure 3**. On the Kromasil column the Clarithromycin peak elutes later at 11.2 minutes. The peak-to-valley ratio between impurity D and clarithromycin on the Kromasil column is 10.0, slightly less than on the Kinetex column.

Conclusion

The results shown above demonstrate that a Kinetex $3.5\,\mu m$ XB-C18, 100×4.6 column can be successfully used to analyze Clarithromycin according to Ph. Eur. Monograph 1651. The peak-to-valley ratio between the API and impurity D for system suitability was 10.8 which is well above the minimum of 3.0 required by the Ph. Eur. monograph method. In comparison to the Kromasil $3.5\,\mu m$ C18, the Kinetex column acquired a faster analysis with slightly improved resolution between clarithromycin and the closely eluting impurity D.

Figure 2. Clarithromycin reference solution D on a Kinetex 3.5 μm XB-C18, 100 x 4.6 (Peak-to-valley ratio is 10.8)


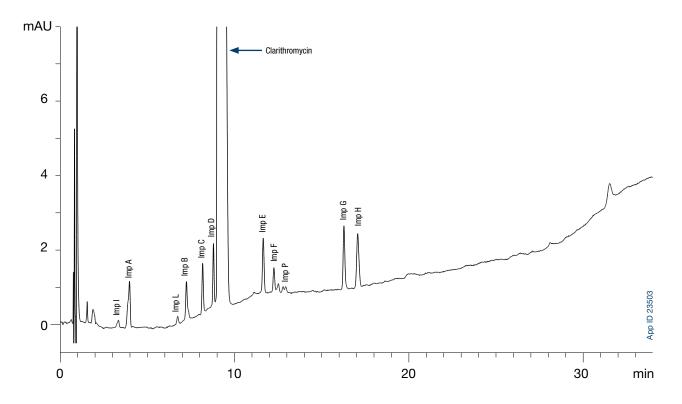




Figure 3. Clarithromycin reference solution D on a Kromasil 3.5 µm C18, 100 x 4.6 (Peak-to-valley ratio is 10.0)

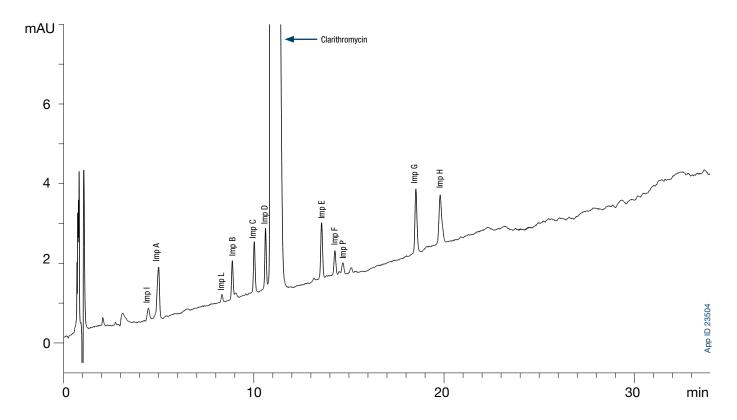

Kinetex 3.5 µm Analytical Columns		SecurityGuard ULTRA Cartridges (mm)
Phases	100 x 4.6 mm	3pk
XB-C18	00F-4744-E0	AJ0-8768
		for 4.6 mm ID

^{*} SecurityGuard ULTRA Cartridges require holder, Part No.: AJ0-9000



If Phenomenex products in this technical note do not provide at least an equivalent separation as compared to other products of the same phase and dimensions, return the product with comparative data within 45 days for a FULL REFUND.

Terms and ConditionsSubject to Phenomenex Standard Terms and Conditions, which may be viewed at http://www.phenomenex.com/TermsAndConditions.

Trademarks

Kinetex is a registered trademark and SecurityGuard is a trademark of Phenomenex. AkzoNobel and Kromasil are registered trademarks of AkzoNobel Pulp and Performance Chemicals AB. Agilent is a registered trademark of Agilent Technologies, Inc.

Disclaimer

Comparative separations may not be representative of all applications.

© 2016 Phenomenex, Inc. All rights reserved.



phenomenex …breaking with tradition[™]

PLICATIONS

Australia

- t: +61 (0)2-9428-6444 f: +61 (0)2-9428-6445 auinfo@phenomenex.com
- Austria
- t: +43 (0)1-319-1301 f: +43 (0)1-319-1300 anfrage@phenomenex.com

Belaium

t: +32 (0)2 503 4015 (French) t: +32 (0)2 511 8666 (Dutch) f: +31 (0)30-2383749 beinfo@phenomenex.com

Canada

t: +1 (800) 543-3681 f: +1 (310) 328-7768 info@phenomenex.com

China

t: +86 (0)20 2282-6668 f: +86 (0)20 2809-8130 chinainfo@phenomenex.com

Denmark

t: +45 4824 8048 f: +45 4810 6265 nordicinfo@phenomenex.com

Finland

t: +358 (0)9 4789 0063 f: +45 4810 6265 nordicinfo@phenomenex.com

France t: +33 (0)1 30 09 21 10 f: +33 (0)1 30 09 21 11 franceinfo@phenomenex.com

Germany t: +49 (0)6021-58830-0 f: +49 (0)6021-58830-11 anfrage@phenomenex.com

t: +91 (0)40-3012 2400 f: +91 (0)40-3012 2411 indiainfo@phenomenex.com

Ireland

t: +353 (0)1 247 5405 f: +44 1625-501796 eireinfo@phenomenex.com

Italy

t: +39 051 6327511 f: +39 051 6327555 italiainfo@phenomenex.com

Luxembourg t: +31 (0)30-2418700 f: +31 (0)30-2383749 nlinfo@phenomenex.com

Mexico t: 001-800-844-5226 f: 001-310-328-7768 tecnicomx@phenomenex.com

The Netherlands t: +31 (0)30-2418700 f: +31 (0)30-2383749 nlinfo@phenomenex.com

New Zealand t: +64 (0)9-4780951 f: +64 (0)9-4780952 nzinfo@phenomenex.com

Norway t: +47 810 02 005

f: +45 4810 6265 nordicinfo@phenomenex.com

Puerto Rico t: +1 (800) 541-HPLC f: +1 (310) 328-7768 info@phenomenex.com

Spain

t: +34 91-413-8613 f: +34 91-413-2290 espinfo@phenomenex.com

Sweden t: +46 (0)8 611 6950 f: +45 4810 6265 nordicinfo@phenomenex.com

United Kingdom

t: +44 (0)1625-501367 f: +44 (0)1625-501796 ukinfo@phenomenex.com

t: +1 (310) 212-0555 f: +1 (310) 328-7768 info@phenomenex.com

All other countries Corporate Office USA t: +1 (310) 212-0555

f: +1 (310) 328-7768 info@phenomenex.com

www.phenomenex.com

Phenomenex products are available worldwide. For the distributor in your country, contact Phenomenex USA, International Department at international@phenomenex.com