

Nitrogen and Phosphorous Pesticides (EPA 507) Mix # 4 on ZB-5ms


Column: Zebron™ ZB-5MS, GC Cap. Column 30 m x 0.25 mm x 0.25 µm, Ea
Phase: 5% Phenyl-Arylene 95% Dimethylpolysiloxane
Dimensions: 30 meters x 0.25 mm x 0.25 µm
Order No: 7HG-G010-11
Oven Profile: 40 °C for 1 min to 100 °C @ 15 °C/min to 225 °C @ 5 °C/min to 320 °C @ 15 °C/min for 8 min
Carrier Gas: Constant Flow Helium, 0.8 mL/min
Injection: On-Column :1 0.5 µL @ 43°C
Detection: Mass Selective (MSD) (335°C)
Analyst Note: Analytes 10 ppm in Acetone



Products used in this application:



17651

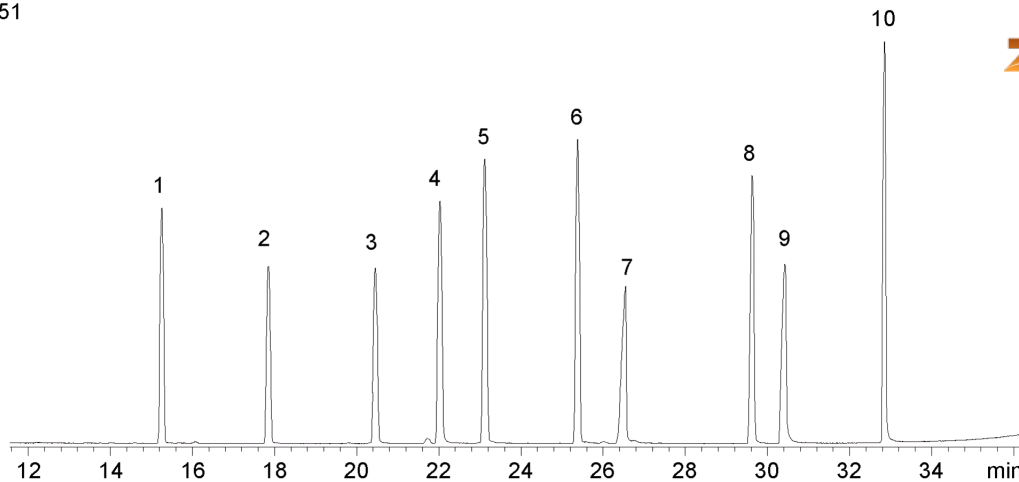

Nitrogen and Phosphorous Pesticides (EPA 507) Mix # 4 on ZB-5ms

ANALYTES:

- 1 Butylate
- 2 Molinate
- 3 Chlorpropham
- 4 Atraton
- 5 Pronamide (Propyzamide)
- 6 Alachlor
- 7 Bromacil
- 8 Tetrachlorvinphos (Tetrachlorovinphos, aka: Stirofos)
- 9 Tricyclazole
- 10 Hexazinone

