

## Nitrogen and Phosphorous Pesticides (EPA 507) Mix # 3 on ZB-5MSi


**Column:** Zebtron™ ZB-5MSi, GC Cap. Column 30 m x 0.25 mm x 0.25 μm, Ea  
**Phase:** 5% Phenyl 95% Dimethylpolysiloxane  
**Dimensions:** 30 meters x 0.25 mm x 0.25 μm  
**Order No:** 7HG-G018-11  
**Oven Profile:** 40 °C for 1 min to 100 °C @ 15 °C/min to 225 °C @ 5 °C/min to 320 °C @ 15 °C/min for 5 min  
**Carrier Gas:** Constant Flow Helium, 0.8 mL/min  
**Injection:** On-Column :1 0.5 μL @ 43°C  
**Detection:** Mass Selective (MSD) (335°C)  
**Analyst Note:** Analytes 10 ppm in Acetone



Products used in this application:



18333

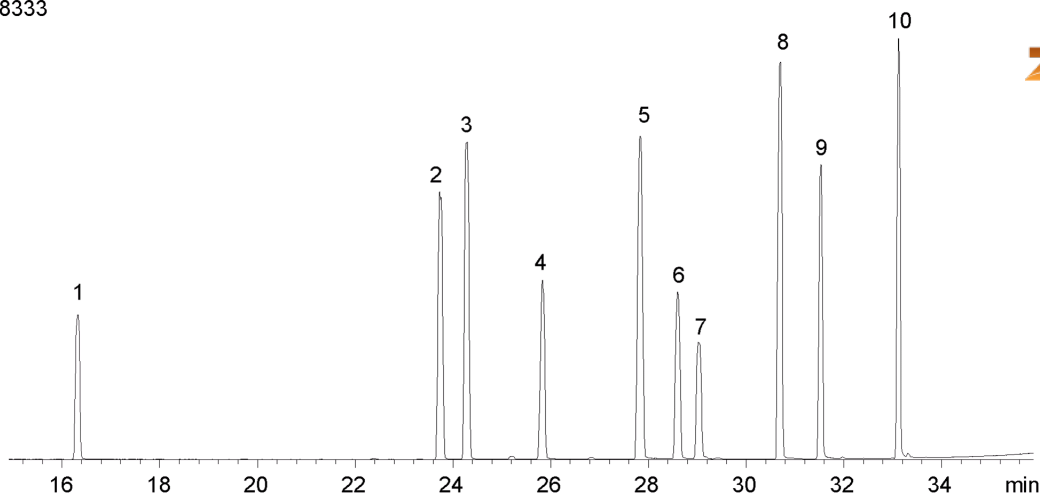

## Nitrogen and Phosphorous Pesticides (EPA 507) Mix # 3 on ZB-5MSi

### ANALYTES:

- 1 Vernolate
- 2 Terbufos (Terbuphos)
- 3 Diazinon
- 4 Metribuzin
- 5 Metolachlor
- 6 MGK-264 (isomer 1)
- 7 MGK-264 (isomer 2)
- 8 Butachlor
- 9 Carboxin
- 10 Norflurazon

